



A Brand Division of American Tooth Industries

1200 Stellar Drive • Oxnard, CA 93033-2404
(805) 487-9868 • Toll Free (800) 235-4639 • FAX: (805) 483-8482
E-mail: info@americantooth.com • Web Site: www.americantooth.com



Justi Cadco Cast Iron Range

General Description:

Our portable range is the solution for high-demand use. The high-grade, long life heating plate is equipment with a variable temperature control. The light will come on ONLY during heat cycle. It will cycle on and off with temperature.

Initial Use:

Heat the appliance without a pressure pot, for approximately 10 minutes, at maximum setting in order for the special coating to seal. Temporary smoking is normal during initial use.

Heating Pot:

This range is only recommended for use with the Justi Pakto® Pressure Pot (070-28-0020).

Directions for Use:

The heating element is equipment with a patented thermostatic system which gives a variable temperature selection. The heating element always heats up with maximum power until it reaches the selected temperature. Thereafter, the thermostat adjusts itself to maintain the temperature; this is indicated by the pilot lamp going on and off.

Maintenance:

The most effective and safe way to clean the range is with a mild detergent and a soft, moist cloth. Do not use any scouring agents.

Warranty:

Six months from the purchase date.



Specifications:

Group/Item#	Description
072-28-0033	Justi Cadco Cast Iron Range
Watts	1500
Amps	12.5
Volts	120, 60 Hz

Note: A short power cord has been provided to reduce the hazards resulting from becoming entangled in or tripping over a long cord. An extension cord may be used if care is properly exercised in its use. The marked electrical rating of the cord should be at least as great as the electrical rating of the range. The longer cord should be arranged so that it will not drape over the counter top or table top where it can be pulled on or tripped over accidentally.

IMPORTANT SAFEGUARDS

To reduce the risk of personal injury or property damage when using electrical appliances, basic safety precautions should always be followed, including the following:

1. Do not touch hot surfaces. Use hot pads or pot holders until unit cools.
2. To protect against electrical shock, do not immerse base, cord or plug into water or other liquid.
3. Unplug from wall outlet when not in use and before cleaning. Allow to cool completely before putting on or taking off parts, and before cleaning or storing the cast iron range.
4. Do not operate with a damaged cord or plug or if unit malfunctions or has been damaged.
5. The use of accessory attachments is not recommended.
6. Do not let cord hang over edge of table or counter or touch hot surfaces.
7. Do not place on or near a hot gas or electric burner or in a heated oven.
8. A fire may occur if unit is covered or touching flammable material including curtains, walls, etc., when in operation.
9. It is recommended that this unit not be moved when it contains a heated pot.
10. Do not clean with scouring pads. They can scratch and ruin the finish of your unit. Pieces can also break off the pad and touch electrical parts, creating a shock hazard.
11. Do not cover the range with metal foils; this may cause overheating of the unit.
12. Do not use this range for other than intended use.

Additional questions, to place your order, or for the Dealer nearest you, please call the Tooth Express Order Desk at 800-628-1437.